PRÁCTICA I.1

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E-mail: pradora@farmacia.ufmg.br

INTRODUCTION

The oil/water partition coefficient (P) is defined for dilute solutions as the concentration ratio of a single species (A) between two phases (organic and aqueous) at equilibrium:


$$P = [A]_{\text{org}} / [A]_{\text{aq}}$$

In recent years, the partition coefficient has become a widely studied and very important physicochemical parameter since its logarithm, usually for the *n*-octanol/water system ($\log P_{\text{o/w}}$) is widely used as a measure of hydrophobicity/lipophilicity. Today, $\log P_{\text{o/w}}$ probably represents the most informative and successful physicochemical property used in medicinal chemistry. It has been extensively employed to explain drug-receptor and drug-biological membrane interactions, and it has been a crucial parameter in many quantitative structure–activity relationships (QSARs) developed for pharmaceutical, environmental, biochemical, or

drug design applications.

The most extensive and useful partition coefficient data were obtained by simply shaking a solute with two immiscible solvents and then analyzing the solute concentration in one of both phases; this is the classical “shake-flask” procedure.

The objective of the practical experiment described here is to determine the partition coefficient of mandelic acid, an antiseptic agent for urinary tract infections, using the shake-flask method. The partition coefficient of mandelic acid described in the literature is 1.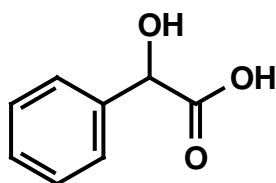

mandelic acid

EXPERIMENTAL

Basic reference: R. T. Smedberg. *J. Chem. Educ.* **71** (3), 269 (1994).

Part 1

- Transfer 15.0 mL of an aqueous solution of mandelic acid (about 1 g/100 mL) to an Erlenmeyer flask, add 20 mL of water, 2 drops of a solution of phenolphthalein indicator (0.1 g in 100 mL of ethanol) and titrate with a standardized solution of sodium hydroxide (about 0.1 mol/L) to a pink color.

Part 2

- Transfer 15.0 mL of the same aqueous solution of mandelic acid used in part 1 (about 1 g/100 mL) to a separatory funnel, add 10 mL of ethyl ether, and shake. Transfer the aqueous layer (bottom) to an Erlenmeyer flask, add 20 mL of water, 2 drops of a solution of phenolphthalein indicator (0.1 g in 100 mL of ethanol) and titrate with a standardized solution of sodium hydroxide (about 0.1 mol/L) to a pink

color.

CALCULATIONS

- Based on the data collected in the part 1, calculate the concentration of mandelic acid in the original solution (mg/mL, g/100 mL, mol/15 mL, mol/10 mL, mol/L).
- Based on the data collected in part 2, calculate the concentration of mandelic acid in the aqueous solution after the extraction with ethyl ether (mg/mL, g/100 mL, mol/15 mL, mol/10 mL, mol/L).
- Based on the data collected in parts 1 and 2, calculate the concentration of mandelic acid in the ether layer (mg/mL, g/100 mL, mol/15 mL, mol/10 mL, mol/L).
- Calculate the partition coefficient of mandelic acid.

OBSERVATION

After the experiment, recover the ethyl ether and the mandelic acid of the organic phase by distillation.

Maria Auxiliadôra Fontes Prado

doraprad@dedalus.lcc.ufmg.br

High standards in safety measures should be maintained in all work carried out in Medicinal Chemistry Laboratories. The handling of electrical instruments, heating elements, glass materials, dissolvents and other inflammable materials does not present a problem if the supervisor's instructions are carefully followed.

This document has been supervised by Prof. Maria Auxiliadôra Fontes Prado) who has informed that no special risk (regarding toxicity, inflammability, explosions),

outside of the standard risks pertaining to a Medicinal Chemistry laboratory exist when performing this exercise.

If your exercise involves any "special" risks, please inform the editor.